

Contec ProChlor SATWIPES System

Meltblown polypropylene wipes with stabilised hypochlorous acid blended with purified water

Contec ProChlor is a revolutionary new sporicide achieving a 100% kill against spores in under 1 min. It is validated for sporicidal transfer disinfection in practice-mimicking conditions, achieving a log 5 kill on syringe paper samples and complete kill on plastic syringe samples, in 1 min.

A patent pending blend of hypochlorous acid in purified water, Contec ProChlor is efficacious against bacteria, fungi, moulds, yeasts, viruses and spores.

Contec ProChlor is provided as a point-of-use SATWIPE system. Each double bag pack contains a canister of 120 x 34gsm meltblown polypropylene wipes and 700ml Contec ProChlor, in a 1L bottle.

Simply add 700ml of ProChlor to the dry roll, 5 mins before use, and use the presaturated wipes within 8 hours.

Contec ProChlor is 0.2 micron filtered, and filled in a Grade A environment. The sterile product is filtered into pre-irradiated components. The dry wipe canister is irradiated at no less than 25 kGy.



FEATURES	BENEFITS
Impregnated wipes at point-of-use	<ul style="list-style-type: none"> Meets requirements of Aseptic Transfer Guidelines
Log 5 kill against <i>Bacillus subtilis</i> in 60 seconds	<ul style="list-style-type: none"> Fast acting so suitable for transfer disinfection
120 polypropylene wipes in a canister	<ul style="list-style-type: none"> Reduces amount of packaging generated Fast and easy to use High quality wipe for transfer disinfection
Filtered to 0.2 microns in a Grade B cleanroom	<ul style="list-style-type: none"> Ensures the product is free from contamination and particulates
Sterile version available	<ul style="list-style-type: none"> Suitable for Grade A and B cleanrooms
Contains no quaternary ammonium or surfactant	<ul style="list-style-type: none"> Very low residue, saving time on residue removal
No hazard classification	<ul style="list-style-type: none"> Good operator acceptability
Double bagged in linear tear packaging	<ul style="list-style-type: none"> Each bag is easy to open even when wearing gloves

For more information or to request a sample,
e-mail infoeu@contecinc.com or phone + 33 (0) 2 97 43 76 98.

Copyright © 2018 Contec, Inc. All rights reserved. INT122 121118

Contec
Vannes Cedex
France
Tel: +33(0) 2 97 43 76 98

Contec Inc
Spartanburg, SC
United States
Tel: +1 (864) 503-8333

Contec Cleanroom Technology (Suzhou) Co. Ltd
Suzhou
China
Tel: +86-512-6274 4050



Product Specification

Contec ProChlor 700ml

200ppm hypochlorous acid in purified water

Ready to use with an 18 month shelf life

0.2 micron filtered in a grade B cleanroom

Aseptic filled under UDAF grade A in a grade B cleanroom

Double bagged in polyethylene linear tear bags

Relevant certificates of analysis and sterility are provided with every batch. Full supporting documentation information is available on request.

SATWIPE Canister 120 wipes

34gsm polyester / cellulose wipes 150 x 200mm

18 month shelf life

Packed in a grade C cleanroom

Gamma irradiated at no less than 25kGy

Efficacy

Contec ProChlor has been shown to have good efficacy against spores using standard laboratory tests. Contec ProChlor SATWIPES have also been tested against the EN 16615 test for presaturated wipes against bacteria, fungi and bacterial spores.

Test	Test Type	Organism	Contact Time	Result	Test	Test Type	Organism	Contact Time	Result **
EN13704	Suspension Test	Bacillus <i>Subtilis</i>	1 min	Log 6	EN16615	Wiping Test	<i>S. aureus</i>	1 min	Log 7.24
EN13704	Suspension Test	Bacillus <i>Cereus</i>	1 min	Log 3.24*	EN16615	Wiping Test	<i>P. aeruginosa</i>	1 min	Log 7.45
EN13704	Suspension Test	Bacillus <i>Pumilis</i>	1 min	Log 3.88*	EN16615	Wiping Test	<i>A. brasiliensis</i>	1 min	Log 6.13
EN13697	Surface Test	Bacillus <i>Subtilis</i>	1 min	Log 6	EN16615	Wiping Test	<i>B. subtilis</i>	1 min	Log 7.11

* Complete kill

** Test fields 2-4 average of 0 cfu.

Product Application

Pour 700ml Contec Sterile ProChlor into the centre of the wipe roll. Close lid and invert canister multiple times for a 20 second period. Wait 5 minutes this will ensure saturation of all wipes. Following the 5-minute saturation time, some liquid will remain in the canister, this is intentional to ensure efficient wetting of the wipes. Once saturated with ProChlor, the wipes are suitable for sporicidal transfer disinfection. Wipes should be used within 8 hours & immediately disposed of. Dispose of remaining fluid and canister according to facility procedures.

Part No.	Description	Size	Packaging
SSWPP02PC	Contec Sterile Meltblown Polypropylene SATWIPES System For use with Contec Sterile ProChlor	150 x 200mm	120 wipes per canister; 10 canisters per case
FSWPP02PC	Contec Meltblown Polypropylene SATWIPES System For use with Contec ProChlor	150 x 200mm	120 wipes per canister; 10 canisters per case

Use biocides safely. Always read the label and product information before use.